UCL010

UCL022

UCL060

HANDLEIDING - MODE D'EMPLOI - MANUAL

UCL010 (754750070)

UCL022 (754750071)

UCL060 (754750073)

Ultrasonne reinigers **Nettoyeurs à ultrasons** **Ultrasonic cleaning machines** **Limpiadores ultrasónicos**

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Warning

Please kindly check this safety instruction when you use this unit, and please keep this manual carefully after reading. Our products series are developing constantly and any technical change, we can adopt without other advise.

General introduction

Ultrasonic cleaning machines, introducing foreign advanced technology and equipment, were developed by technologists through many years of researches.

The ultrasonic cleaning machine made by our factory is widely used in fields like hospitals, pharmacies, schools, scientific researches, petroleum, chemical industry, light industry, metallurgy, machines, transportations, national defences etc. Currently it's an ideal, efficient and accurate cleaning equipment to clean thoroughly the instrument, meter, electronic device, circuit board, magnetic material, galvanized parts, hardware, optical lens, audio-frequency magnetic head, spinneret, latex mold, medical apparatus, glassware, bijouterie, parts of a clock and watch, fine metal, bearing, nozzle, oil pump and parts in manufacturing machine and some shape-complicated parts, especially the deep hole, blind hole and concave or convex surfaces.

Working principle

The ultra-audio-frequency electric signal produced by the ultrasonic generator in a ultrasonic cleaning machine is converted to the mechanical vibration with the same frequency through the piezoelectric inverse-transforming effect of the transducer, and then it will radiate forward intensively in the cleaning solution in the form of ultra-audio-frequency longitudinal wave (The ultra-audio frequency longitudinal wave spreads alternatively between zones of negative pressure and positive pressure), producing ten thousands of small bubbles to explode in turn. It's these small bubbles which form and grow in the zone of negative pressure, and explode instantaneously in the zone of positive pressure, to strike with high pressure the surface, gap, and small local of a cavity to be washed, leading to peeling off of fouling quickly, thus the good cleaning and purifying effect can be reached.

Safety instructions

1. Before using, please read the manual carefully.
2. Before using, make sure the ground connection should be OK.
3. No fire can be permitted in the circumstance.
4. When using the liquid which is easy to be fired, please take care of it.
5. Please don't let the temperature of liquid over 80°C, generally speaking, normal and middle temperature is enough.
6. If no liquid in the tank, don't start the machine and heating, avoid the machine damaged under no-loading condition.
7. You should use good-quality, big power plug, don't use bad quality plug, don't use this together with air conditioner and electric heater, avoid the current too big.
8. Only personality can be permitted to use it.
9. Don't use big power unit to be working together with this machine, don't use this machine into the water to working.

Technical specifications

Model	UCL 010	UCL 022	UCL 060
Capacity	10 l	22 l	60 l
Ultrasonic power	240 W	500 W	1000 W
Heating power	285 W	475 W	1500 W
Total power	525 W	975 W	2500 W
Inner chamber size	300 x 240 x 150 mm	500 x 300 x 150 mm	500 x 400 x 300 mm
Frequency	40 kHz	40 kHz	28 kHz
Adjustable temperature	~ 80°C	~ 80°C	~ 80°C
Adjustable time	1 ~ 99 min	1 ~ 99 min	1 ~ 99 min
Drainage	Yes	Yes	Yes
Basket	Yes	Yes	Yes
Lower-noise cover	Yes	Yes	Yes

Packing list

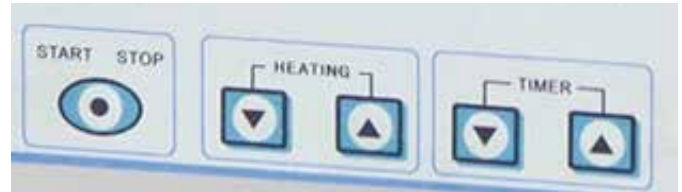
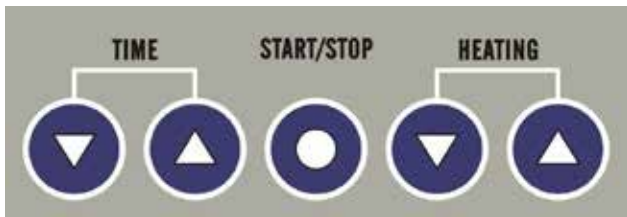
- 1 Cleaning machine
- 1 Operation manual
- 1 Q.C. Pass
- 1 Basket
- 1 Power line

Note: No drainage and drainage pipe for under 5l

Note: The technology and specifications can be altered by us without notice.

Instructions for use

1. Check the grounding wire of the ultrasonic cleaning machine and power line is in good condition (electricity leaking exists or not).
2. The cleaning solution can be put into the cleaning slot. When cleaning, level of the solution can't be below one third of the height of the cleaning slot, above two thirds of the cleaning slot. You should add different detergents according to different cleaning requirements, thus to increase cleaning efficiency. All detergents should not fret the inferior slot and the body of the cleaning machine.
3. Insert the plug of the ultrasonic cleaning machine into the power socket of 220-240VAC - 100-120VAC. Rotate the switch of the timing controller and the indicator will be on. When the surface of the solution fluctuates like a spider web with noise, it means that the cleaning machine is working.
4. According to the cleaning requirement, set the cleaning temperature.
5. Put things to be washed into the metal frame, and set the cleaning time according to the real situation of fouling accumulation. Normally the time can be set between 3-30 minutes, and the time can be prolonged for things which are especially difficult to be cleaned.
6. For heavy articles, suspend it through a hanger in the cleaning solution to increase the cleaning effect, heighten the cleaning quality.
7. Take the article to be washed out; rewash or rinse with the ultrasonic machine.
8. If you don't use the machine for a long time, you should drain the cleaning solution in the slot out, and clean the body of the machine.

Temperature-controlling timing on the LED

1. Set the ultrasonic cleaning time (99 minutes).
2. Press "increase" or "deduce" to set needed cleaning time.
3. Set the temperature (~80°C).
4. Press "increase" or "deduce" to set the temperature of the needed cleaning solution.
5. Press "On/off" key to start the machine to work.

Service and maintenance

This machine is very valuable, so a special man should be assigned to be responsible for operation, service and maintenance. Thorough checking to the machine should be done periodically.

1. Check the looseness, overheating, dampness and the contact of every connection of the power and the output line.
2. Check the mesh enclosure of the power cooling fan, and please remove the dust on it.
3. Check the leakage of the sealing of the supporting plate.
4. Please contact the supplier if the ultrasonic oscillator falls off or the power tube is burned.

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